

APPLICATIONS

Determination of Perfluoroalkyl Substances (PFAS) in Sediments by QuEChERS Extraction and HPLC-MS/MS

Syljohn Estil¹, Eric Nelson¹ and Scott Krepich²

¹Sanitation Districts of Los Angeles County-San Jose Creek Water Quality Laboratory, Whittier CA

²Phenomenex, Inc., 411 Madrid Avenue., Torrance, CA 90501



Scott Krepich
Senior Field Application
Scientist

Scott enjoys surfing and eating. He is crazy about chromatography, because his mom is really into CSI and thinks that is what he does.

Introduction

Perfluoroalkyl substances (PFAS) are a class of highly stable synthetic organic compounds used in a wide variety of industrial and commercial applications including surface treatment for textiles, packaging materials, and non-stick cookware. PFAS are characterized by a hydrophobic fully fluorinated alkyl chain and a hydrophilic functional group. They are persistent in the environment due to the exceptional stability of the C-F bond. Once released in the aquatic environment, these chemicals will partition between the water phase and the sediment. To better understand the fate and transport of these compounds, it is important to measure both the solid and liquid environmental fractions.

There are several methods available for the extraction and analysis of PFAS in aqueous samples. However, very few procedures are available for extracting these compounds in solid matrices such as sediments. Typical methods used are mechanical shaker and ultrasonic-assisted Solid-Liquid Extractions (SLE)^{6,7,8}. The extracts are then subjected to additional cleanup steps, usually by solid phase extraction. These are generally solvent-intensive and time-consuming processes. In 2003, an extraction procedure called QuEChERS (Quick-Easy-Cheap-Effective-Rugged-and-Safe) was introduced. It was originally developed to extract pesticide residues in food matrices but has since found applications in the field of environmental analytical chemistry.

Our laboratory (LACSD) previously developed and validated a QuEChERS sediment extraction procedure for emerging contaminants including: pharmaceutical and personal care products, steroids, alkylphenol ethoxylates, and pyrethroid pesticides^{2,3,4}. We have successfully applied the same extraction method to determine perfluoroalkyl substances in marine and freshwater sediments.

Materials and Methods

Reagents/Chemicals

- QuEChERS Extraction – In a 50 mL plastic centrifuge tube combine 2.0 g of Anhydrous Magnesium Sulfate, and 1.5 g Sodium Acetate or use approximately 3.5 g of AOAC2007.01 roQ™ extraction packet (part no. AH0-9043)
- QuEChERS dSPE Clean-Up – roQ 15 mL dSPE Kit (part no. KS0-8926)

Sample Preparation

QuEChERS Extraction Protocol

1. Weigh 2.0 g of dried sediment into a polypropylene container and spike with isotopically-labeled internal standards. PPCPs, Steroids, and Pyrethroids can be extracted concurrently with this method by adding the appropriate internal standard and spiking solutions to the samples and QCs^{2,3,4}.
2. Add 10mL deionized water and vortex. Add 10mL acidified acetonitrile (1% acetic acid) to the slurry and vortex.
3. Add the extraction salts (1.5 g Sodium Acetate and 2 g MgSO₄) to the sample and vortex for 1 minute.
4. Centrifuge the samples for 5 minutes at 4000 rpm.
5. Place the samples in a rack and freeze at -20° for 30-60 minutes. This freezing step allows for easier extraction of the supernatant.
6. Transfer 8-9 mL of the acetonitrile supernatant into a roQ QuEChERS PSA/C18 dSPE clean-up tube (Part no. KS0-8926) and vortex for one minute.
7. Centrifuge the dSPE tubes for 10 minutes at 3000 rpm.
8. Place an aliquot of the extract in a HPLC vial and dilute 1:1 with deionized water. The sample is now ready for analysis.

HPLC-MS/MS Conditions

Column:	Gemini® 3 µm C18
Dimensions:	100 x 3 mm
Part No.:	00D-4439-Y0
Inline Filter:	Phenomenex Krudkatcher™ Ultra
Delay Column:	Luna® 5 µm C18 (2) 30 x 2.0 mm
Part No.:	00A-4252-B0
Mobile Phase:	A: 20 mM Ammonium acetate in water B: Methanol
Gradient:	Time (min) % B
	0.0 10
	1.5 65
	8.0 95
	8.1 99
	12.0 99
	12.5 10
Injection:	90 µl
Flow Rate:	0.6 mL/min
Temperature:	40 °C
Detector:	SCIEX 5500 QTRAP®
Detection:	MS/MS ESI Negative (sMRM)
Analytes:	1. PFBA 10. PFOS
	2. PFPeA 11. PFNA
	3. PFBS 12. PFOSA
	4. PFHxA 13. PFNS
	5. PFPS 14. PFDA
	6. PFHxS 15. PFDS
	7. PFHpA 16. PFUdA
	8. PFHpS 17. PFDaA
	9. PFOA 18. PFTrDA
	19. PFTeDA



Mass Spectrometer Parameters

Table 1.
MRM Transitions and Compound Dependent Parameters

Compound Name	Q1	Q3	DP	CE
Perfluorobutanoic acid (PFBA)	213	169	-71	-14
Perfluoropentanoic acid (PFPeA)	263	219	-71	-12
Perfluorohexanoic acid (PFHxA)	313	269	-60	-14
Perfluoroheptanoic acid (PFHpA)	363	319	-62	-15
Perfluorooctanoic acid (PFOA)	413	369	-91	-12
Perfluorononanoic acid (PFNA)	463	419	-79	-15
Perfluorodecanoic acid (PFDA)	513	469	-83	-17
Perfluoroundecanoic acid (PFUdA)	563	519	-60	-17
Perfluorododecanoic acid (PFDoA)	613	569	-50	-21
Perfluorotridecanoic acid (PFTrDA)	663	619	-49	-18
Perfluorotetradecanoic acid (PFTeDA)	713	669	-63	-20
Perfluorobutanesulfonate (PFBS)	299	80	-94	-70
Perfluoropentanesulfonate (PFPeS)	349	80	-96	-66
Perfluorohexanesulfonate (PFHxS)	399	80	-92	-75
Perfluoroheptanesulfonate (PFHpS)	449	80	-75	-84
Perfluorooctanesulfonate (PFOS)	499	80	-78	-96
Perfluorooctanesulfonamide (PFOSA)	498	78	-60	-84
Perfluorononanesulfonate (PFNS)	549	80	-87	-100
Perfluorodecanesulfonate (PFDS)	599	80	-55	-100
Perfluoro-n- ¹³ C4-butanoic acid (M4PFBA)	217	171.9	-71	-13
Perfluoro-n- ¹³ C5-pentanoic acid (M5PFPeA)	268	222.7	-71	-12
Perfluoro-n- ¹³ C5-hexanoic acid (M5PFHxA)	318	272.9	-60	-13
Perfluoro-n- ¹³ C4-heptanoic acid (M4PFHpA)	367	321.8	-62	-14
Perfluoro-n- ¹³ C8-octanoic acid (M8PFOA)	421	376	-91	-12
Perfluoro-n- ¹³ C9-nonanoic acid (M9PFNA)	472	427	-79	-17
Perfluoro-n- ¹³ C6-decanoic acid (M6PFDA)	519	474	-83	-21
Perfluoro-n- ¹³ C7-undecanoic acid (M7PFUdA)	570	525	-60	-17
Perfluoro-n- ¹³ C2-dodecanoic acid (M2PFDoA)	615	570	-50	-24
Perfluoro-n- ¹³ C2-tetradecanoic acid (M2PFTeDA)	715	670	-63	-25
Perfluoro- ¹³ C3-butanesulfonate (M3PFBS)	302	80	-94	-55
Perfluoro- ¹³ C3-hexanesulfonate (M3PFHxS)	402	80	-92	-85
Perfluoro- ¹³ C8-octanesulfonate (M8PFOS)	507	80	-78	-100

Note: DP = Declustering Potential

CE = Collision Energy


Table 2.
MS Source Parameters

Source Parameters	Settings
Temperature	400 °C
Gas 1	50
Gas 2	50
Curtain Gas	35
Ionization Energies	-4500 V
Collision Gas	High

Table 3.
Method Performance Data for Sediments Spiked at 1 ng/g of the Target Analytes (n=4)

Compound	Average % Recovery	% RSD
PFBA	91.7	0.76
PFPeA	86.3	6
PFHxA	89.4	1.2
PFHpA	93.1	2.9
PFOA	98.3	1.5
PFNA	93	1.6
PFDA	87.7	4.5
PFUdA	92.3	2.1
PFDoA	92.5	4.1
PFTrDA	88.2	2.1
PFTeDA	87.6	2.1
PFBS	86.3	2.1
PFPeS	96.2	3.2
PFHxS	81.3	5
PFHpS	92.3	2.6
PFOS	92.1	2.6
PFOSA	104.5	6.3
PFNS	89.8	6.8
PFDS	87.3	6.7

Figure 1
 Extracted ion chromatogram of sediments spiked with 1.0 ng/g of the target analytes

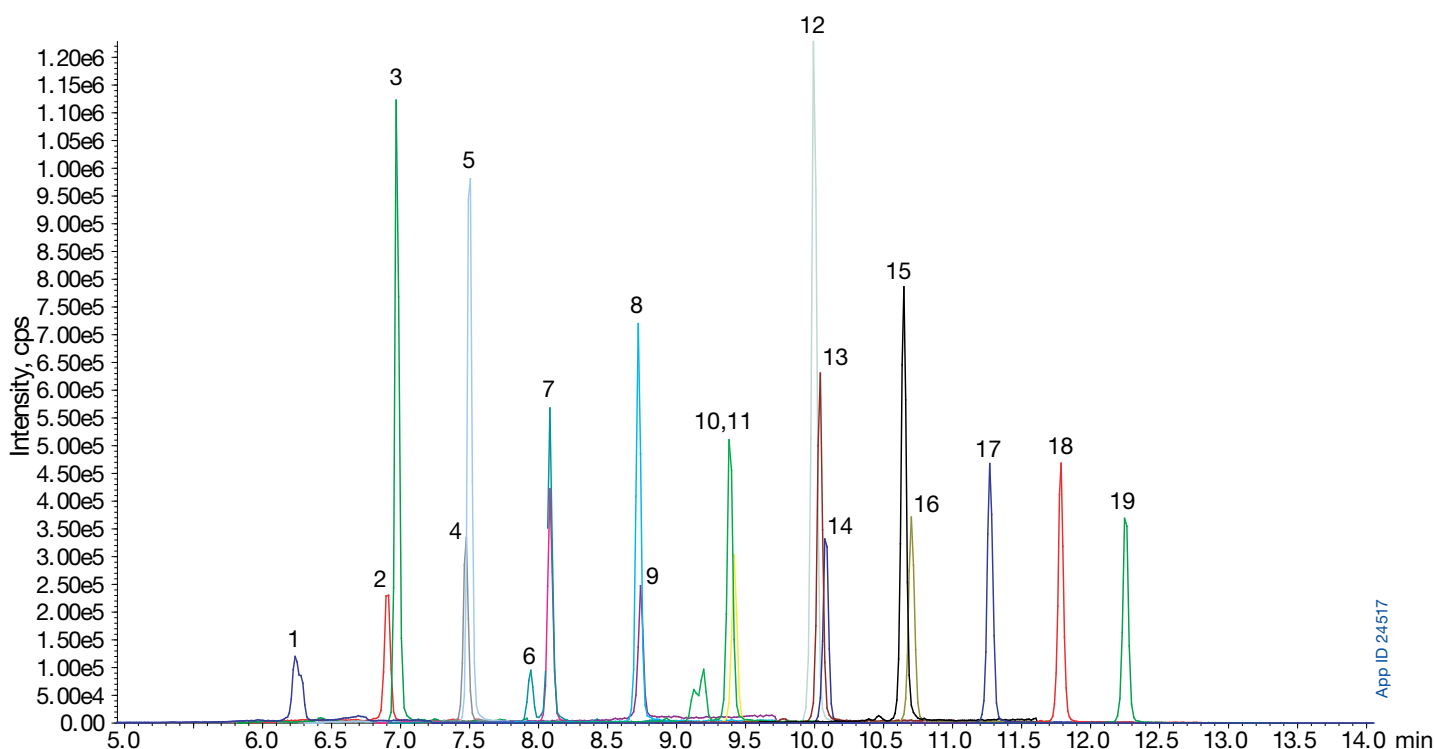

Results and Discussion

QuEChERS is a vortex-assisted solid-liquid extraction procedure that uses acetonitrile, salts, and buffering agents for extraction, phase-separation, and pH adjustment respectively. Extracts are subsequently transferred to a dispersive solid phase extraction (dSPE) tube containing a drying agent ($MgSO_4$) and SPE sorbents such as C18 or PSA for sample cleanup.

The modified QuEChERS method presented here is a simple, efficient, and cost-effective method for determining PFAS levels in sediments. Accuracy and precision were assessed using four replicates of sediments spiked with the target analytes. Average % recoveries are all within the 80-120% range and % RSDs for all analytes are below 10% (**Table 3**). Reporting limits were set at 0.05 ng/g dry weight based on a 2.0 g initial sample weight.

Conclusion

The sediment extraction using the modified QuEChERS method is a fast, effective, and efficient way of extracting 19 perfluoroalkyl substances from marine and river sediment matrices. The procedure significantly minimizes sample preparation time, solvent consumption, and overall cost of analysis. Only minor modifications were made to the well-established QuEChERS method to accommodate the target analytes. Excellent extraction recoveries and precision were achieved and method reporting limits are in the low ng/g range.

References

- Berlitz-Barbier, A., et al., Multi-residue analysis of emerging pollutants in sediment using QuEChERS-based extraction followed by LC-MS/MS analysis. *Analytical and Bioanalytical Chemistry* (2014) 406:1259-1266
- Estil S, et al., A. Rapid Extraction and Analysis of PPCPs from Sediments by QuEChERS and LC/MS/MS (2016) Phenomenex Application Note: TN-0099
- Estil S, et al., A. Rapid Extraction and Analysis of Steroids and Pyrethroids from Sediments by QuEChERS and LC/MS/MS (2016) Phenomenex Application Note: TN-0096 & TN-0098
- Estil S, et al., A. Rapid Extraction and Analysis of Pyrethroids from Sediments by QuEChERS and LC/MS/MS (2016) Phenomenex Application Note: TN-0098
- Roberts S, et al. Quantitation of PFAS in Water Samples using LC/MS/MS: Large Volume Direct Injection and Solid Phase Extraction (2016) SCIEX Application Note Publication Number: RUO-MKT-02-4707-A
- Jahnke A, et al. Trace analysis of per- and polyfluorinated alkyl substances in various matrices – How do current methods perform? *Journal of Chromatography A* (2009) 1216: 410-421
- Ahrens et. al. Partitioning behavior of per- and polyfluoroalkyl compounds between pore water and sediment in two sediment cores from Tokyo Bay, Japan. *Environmental Science and Technology* (2009) 43: 6969-6975
- Long et. al. Quantification of Pharmaceuticals, Personal Care Products, and Perfluoroalkyl Substances in the Marine Sediments of Puget Sound, Washington, USA. *Environmental Toxicology and Chemistry* (2013) 32: 1701-1710

Acknowledgements

Special thanks Syljohn Estil and to the Sanitation Districts of Los Angeles County – San Jose Creek Water Quality Laboratory for contributing this method.



APPLICATIONS

Gemini[®] HPLC Columns Ordering Information

3 μ m Microbore, Minibore and MidBore [™] Columns (mm)										SecurityGuard [™] Cartridges (mm)	
Phases	50 x 1.0	20 x 2.0	30 x 2.0	50 x 2.0	100 x 2.0	150 x 2.0	50 x 3.0	100 x 3.0	150 x 3.0	4 x 2.0*	
Gemini [®] C18	00B-4439-A0	00M-4439-B0	00A-4439-B0	00B-4439-B0	00D-4439-B0	00F-4439-B0	00B-4439-Y0	00D-4439-Y0	00F-4439-Y0	AJ0-7596	

for ID: 2.0-3.0 mm

Luna[®] HPLC Columns Ordering Information

5 μ m Microbore and Minibore Columns (mm)								SecurityGuard [™] Cartridges (mm)	
Phases	50 x 1.0	150 x 1.0	250 x 1.0	30 x 2.0	50 x 2.0	150 x 2.0	250 x 2.0	4 x 2.0*	
Luna [®] C18(2)	00B-4252-A0	00F-4252-A0	00G-4252-A0	00A-4252-B0	00B-4252-B0	00F-4252-B0	00G-4252-B0	AJ0-4286	

for ID: 2.0-3.0 mm

Australia

t: +61 (0)2-9428-6444
f: +61 (0)2-9428-6445
auinfo@phenomenex.com

Austria

t: +43 (0)1-319-1301
f: +43 (0)1-319-1300
anfrage@phenomenex.com

Belgium

t: +32 (0)2 503 4015 (French)
t: +32 (0)2 511 8666 (Dutch)
f: +31 (0)30-2383749
beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681
f: +1 (310) 328-7768
info@phenomenex.com

China

t: +86 400-606-8099
f: +86 (0)22 2532-1033
phen@agela.com

Denmark

t: +45 4824 8048
f: +45 4810 6265
nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063
f: +45 4810 6265
nordicinfo@phenomenex.com

France

t: +33 (0)1 30 09 21 10
f: +33 (0)1 30 09 21 11
franceinfo@phenomenex.com

Germany

t: +49 (0)6021-58830-0
f: +49 (0)6021-58830-11
anfrage@phenomenex.com

India

t: +91 (0)40-3012 2400
f: +91 (0)40-3012 2411
indiainfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405
f: +44 1625-501796
eireinfo@phenomenex.com

Italy

t: +39 051 6327511
f: +39 051 6327555
italiainfo@phenomenex.com

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department at international@phenomenex.com

Luxembourg

t: +31 (0)30-2418700
f: +31 (0)30-2383749
nlinfo@phenomenex.com

Mexico

t: 01-800-844-5226
f: 001-310-328-7768
tecnicomx@phenomenex.com

The Netherlands

t: +31 (0)30-2418700
f: +31 (0)30-2383749
nlinfo@phenomenex.com

New Zealand

t: +64 (0)9-4780951
f: +64 (0)9-4780952
nzinfo@phenomenex.com

Norway

t: +47 810 02 005
f: +45 4810 6265
nordicinfo@phenomenex.com

Puerto Rico

t: +1 (800) 541-HPLC
f: +1 (310) 328-7768
info@phenomenex.com

Spain

t: +34 91-413-8613
f: +34 91-413-2290
esinfo@phenomenex.com

Sweden

t: +46 (0)8 611 6950
f: +45 4810 6265
nordicinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367
f: +44 (0)1625-501796
ukinfo@phenomenex.com

USA

t: +1 (310) 212-0555
f: +1 (310) 328-7768
info@phenomenex.com

All other countries Corporate Office USA

t: +1 (310) 212-0555
f: +1 (310) 328-7768
info@phenomenex.com

roQ[™] Extraction Kits

Extraction Kits contain fifty easy-pour salt packets and fifty 50 mL stand-alone centrifuge tubes

Description	Unit	Part No.
EN 15662 Method Extraction Kits		
4.0 g MgSO ₄ , 1.0 g NaCl, 1.0 g SCTD, 0.5 g SCDS	50/pk	KS0-8909*
AOAC 2007.01 Method Extraction Kits		
6.0 g MgSO ₄ , 1.5 g NaOAc	50/pk	KS0-8911*
Original Non-Buffered Method Extraction Kits		
4.0 g MgSO ₄ , 1.0 g NaCl	50/pk	KS0-8910
6.0 g MgSO ₄ , 1.5 g NaCl	50/pk	KS0-8912

*AOAC and EN Extraction Kits also available in traditional non-collared 50 mL centrifuge tubes, Part No.: KS0-8911-NC and KS0-8909-NC

roQ dSPE Kits

dSPE Kits contain pre-weighed sorbents/salts inside 2 mL or 15 mL centrifuge tubes

Description	Unit	Part No.
2 mL dSPE Kits		
150 mg MgSO ₄ , 25 mg PSA, 25 mg C18-E	100/pk	KS0-8913
150 mg MgSO ₄ , 25 mg PSA, 2.5 mg GCB	100/pk	KS0-8914
150 mg, MgSO ₄ , 25 mg PSA, 7.5 mg GCB	100/pk	KS0-8915
150 mg MgSO ₄ , 25 mg PSA	100/pk	KS0-8916
150 mg MgSO ₄ , 50 mg PSA, 50 mg C18-E, 50 mg GCB	100/pk	KS0-8917
150 mg MgSO ₄ , 50 mg PSA, 50 mg C18-E	100/pk	KS0-8918
150 mg MgSO ₄ , 50 mg PSA, 50 mg GCB	100/pk	KS0-8919
150 mg MgSO ₄ , 50 mg PSA	100/pk	KS0-8920
15 mL dSPE Kits		
900 mg MgSO ₄ , 150 mg PSA, 150 mg C18-E	50/pk	KS0-8921
900 mg MgSO ₄ , 150 mg PSA, 15 mg GCB	50/pk	KS0-8922
900 mg MgSO ₄ , 150 mg PSA, 45 mg GCB	50/pk	KS0-8923
900 mg MgSO ₄ , 150 mg PSA	50/pk	KS0-8924
1200 mg MgSO ₄ , 400 mg PSA, 400 mg C18-E, 400 mg GCB	50/pk	KS0-8925
1200 mg MgSO ₄ , 400 mg PSA, 400 mg C18-E	50/pk	KS0-8926
1200 mg MgSO ₄ , 400 mg PSA, 400 mg GCB	50/pk	KS0-8927
1200 mg MgSO ₄ , 400 mg PSA	50/pk	KS0-8928

Bulk roQ QuEChERS Sorbents

Phases	10 g	100 g
C18-E	—	04G-4348
GCB (Graphitized Carbon Black)	04D-4615	04G-4615
PSA	—	04G-4610



If Phenomenex products in this technical note do not provide at least an equivalent separation as compared to other products of the same phase and dimensions, return the product with comparative data within 45 days for a FULL REFUND.

Terms and Conditions

Subject to Phenomenex Standard Terms and Conditions which may be viewed at www.phenomenex.com/TermsAndConditions.

Trademarks

Luna and Gemini are registered trademarks, KrudKatcher, roQ, MidBore and SecurityGuard are trademarks of Phenomenex. QTRAP is a registered trademark of AB SCIEX Pte. Ltd. AB SCIEX[™] is being used under license. SecurityGuard is patented.

© 2017 Phenomenex, Inc. All rights reserved.